

ward's science+ DataHub

The Power of Twelve
The Convenience of One

Getting to Know the DataHub



Chromebook Version



wardsci.com























How to....

Set up the Datahub from the Datahub Menu	1
Pair your Datahub with a Chromebook Laptop via Bluetooth.....	1
Connect the Datahub to a Chromebook Laptop Using a Bluetooth Connection.....	2
Set up the Datahub from the Datahub Menu Using a Bluetooth Connection.....	3
Download Your Data	3
Clear the Datahub Memory	3
Use the Datahub Software (Chromebook Laptop)	4
The Menu/Tool Bar	4
Open Data Files	4
Save	4
Export	4
Statistics	4
Set up	4
Run/Stop	4
Download	4
Mathematical Functions	4
Activities	5
Other Software Features.....	5
Zoom	4
Marker.....	5
Annotate	5
Display Options	5
Use the Datahub Sensors.....	6
Accelerometer.....	6
Air Pressure	6
Ambient Temperature Sensor.....	6
Barometric Pressure Sensor.....	6
Colorimeter	6
Conductivity Sensor	7
Current Sensor	7
Dissolved O2 Sensor.....	7
Distance/Motion Sensor	7
External Temperature Sensor	7
GPS (Global Positioning Satellite) Sensor.....	8
Heart Rate Sensor	8

Humidity Sensor	8
Infrared Sensor.....	8
Light Sensor	8
Microphone.....	9
pH Sensor	9
Sound Sensor	9
Thermocouple	9
Turbidity Sensor	10
UV Sensor	10
Voltage	10
Available External Sensors	10
Dymo Wireless Force & Accelerometer	10
Voltage Sensor	11
CO2 Sensor	11
Heart Rate/Pulse	11
Respiration	11
Current	11
Magnetic Field.....	12

How To...

Set up the Datahub from the Datahub Menu

1. **Turn on** the Datahub by pressing the On/Off key. 
2. **Rotate the Datahub plastic ring** to expose the sensors. Connect any external sensors that you might be using (pH, Dissolved O2, Conductivity, Thermocouple, External Temperature, Force/Accelerometer, Heart Rate, Current, Voltage or Magnetic Field).
3. **Select Sensors** - Press Scroll key  to open the sensor menu. Press the Select key  to open the Set Up menu  and then choose the Set Sensor icon  using the Select  key again. Select the sensors that you wish to use by pushing the buttons on the perimeter of the Datahub. Some buttons control more than one sensor. Push the button multiple times to scroll through the available options. When the sensor(s) that you want is/are highlighted push the Escape key  to return to the set up menu. *Note: Some buttons can only activate one sensor at a time because of the external connections needed by the sensors.*
4. **Set Sampling Rate** - Press the Escape key  to return to the previous menu and use the scroll key  to move to the Sampling Rate icon  and push the Select key . Use the Scroll key  to move through the options. Use the Select key  to choose the rate at which you want to select samples. Press the Select key to confirm your choice. Options include manual, 1/second, 10/second, 100/second, 1000/second and 25000/second. The rate of collection by some sensors is limited and so you may not see all of the options for every sensor.  *Note if you choose to sample manually, skip Step #5.*
5. **Set the Number of Samples** - Press the Escape key  to leave the previous menu and use the scroll key  to move to the Number of Samples icon  and push the Select key.  Use the Scroll key  to choose the number of samples you wish to collect. Options include 10, 100, 1000, 10000. Press the Select key  to confirm your choice. Press the Escape key  three times (3x) to exit the set up menu
6. **Begin Collecting Data** - When you are ready, press the Select key  to begin the data collection.

Pairing your Datahub with a Chromebook Laptop via Bluetooth. The Chromebook version is called Globilab.


1. Download and install the Globilab app from the Google Play Store at:
<https://chrome.google.com/webstore/detail/Datahub/caelfgcgafddnfhocinleaceplgdddpb>
2. Turn on your Datahub.


3. Click on the Chromebook bottom right menu bar




The below dialog box opens:


How To...

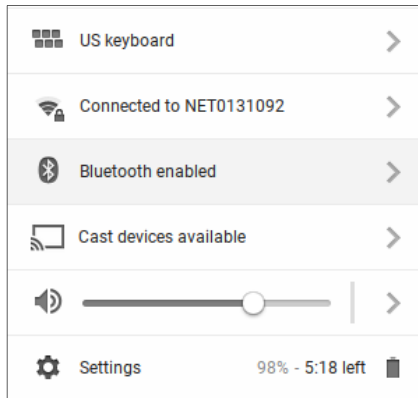

4. Click on the Bluetooth icon. Make sure that Bluetooth is enabled on your Chromebook. Click on **Manage Devices...** then click **Add Device**. The Chromebook will start searching for new Datahub devices.
5. The last 4 digits of your Datahub's serial number will appear on the Chromebook screen as: **Datahub-XXXX** - where the XXXX represent the Datahub last four S/N digits.
6. From the Chromebook's list of devices, select your Datahub. Make sure that the four digit code on the back of the Datahub matches the one you have selected on your Chromebook. Click **Connect**.
7. Press and hold the Scroll key  until you hear a long beep and the Datahub screen reads BT Pairing.
8. On the Chromebook use the PIN code **1234** to pair your Datahub to your Chromebook.
9. Your Datahub should be added to the Chromebook list of paired devices. On some Chromebooks you might see an error message. You can ignore it, as long as you see the Datahub on the Paired devices list.

Connect the Datahub to a Chromebook Laptop Using a Bluetooth Connection


1. Open the Datahub software and turn on the Datahub by pressing the On/Off key. 
2. Make sure that your Datahub has been paired with your computer. (see above) You will only have to pair the Datahub once, the first time you use your Datahub with the Chromebook.
3. From the Datahub menu select "Connect Datahub".

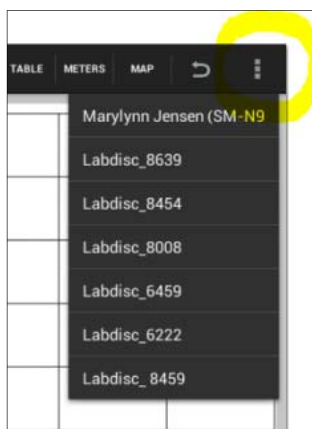

How To...


4. Select the Datahub that you want to connect to from the pop up menu. Make sure the 4 digit code matches the one on the back of the Datahub.


Set up the Datahub from the Datahub Menu Using a Bluetooth Connection

1. Press the “Logger Setup” button  to get the Logger Setup window. Different sensors will show for different Datahub models, depending on what sensors are available. If you are using a Globisens brand external sensor attached to the universal adapter port, it will appear in the set-up menu after you have attached it.
2. Click the check box next to the sensor name to select it. Click again to deselect it.
3. Slide your mouse pointer up and down next to Rate to select the sample frequency.
4. Slide your mouse pointer up and down next to Samples to select the number of samples you wish to collect.
5. Click back on the main screen to exit the Setup window.

Logger Setup	
<input type="checkbox"/> UV	1/Min
<input type="checkbox"/> pH	Rate 1/Sec
<input type="checkbox"/> Barometer	10/Sec
<input type="checkbox"/> Sound	
<input type="checkbox"/> IR Temperature	
<input type="checkbox"/> Humidity	1000
<input type="checkbox"/> GPS	Samples 10000
<input checked="" type="checkbox"/> Amb. Temperature	10
<input checked="" type="checkbox"/> Ext. Temperature	
<input type="checkbox"/> Colorimeter	
<input type="checkbox"/> Turbidity	
<input type="checkbox"/> DO ₂ Dissolved oxygen	
<input type="checkbox"/> Ext. Analog sensor	

Download Your Data

1. Make sure your Datahub is connected to your Chromebook and that it is not running/collecting data.
2. Click the download button  to download a specific data collection set into the Datahub software. This will allow you to select from a list of prior data collection sets.

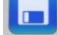Select the data set that you wish to download. Save the data collection by Clicking the save button.  Be sure to clearly label your data collection so that you will not confuse it with the other data collections.

Clear the Datahub Memory













1. Make sure your Datahub is connected to your Chromebook and that it is not running/collecting data.
2. From the Datahub menu select “Clean Datahub”.

METERS	MAP	↶	⋮
Clean Labdisc			
Disconnect Labdisc_4176			
Connect Labdisc			
Settings			


How To...

Use the Datahub Software (Chromebook Laptop)

The Menu/Tool Bar

	Open Data Files – open graphs showing experiment data previously collected and saved. This includes both sample data files provided with the Datahub software and data that you have collected.
	Save – save data from a current or downloaded experiment. Be sure to label your files clearly for future use.
	Export – export the data in chart form to spreadsheet.
	Statistics – quickly see the maximum and minimum values for a specific data collection as well as the average reading, the standard deviation and the number of samples and sampling rate.
	Set up – select which sensors and set up the desired collection rate and collection size from the Chromebook rather than from the Datahub itself. In order to use the set up button, the Datahub must be connected to the Chromebook and must not be running an experiment. Keep in mind that different sensors have different maximum sampling rates. Datahub will limit the collection rate to that of the slowest sensor.
	Run/Stop – begin or stop data collection. Like the set up button, this button will only work if the Datahub is connected to the Chromebook.
	Download – select and download data from any experiment remaining on the Datahub. The Datahub's memory can save up to 127 recordings and will show date/time, number of samples, rate of collection and which sensors were used.
	Zoom – enlarge a section of the graph. <ul style="list-style-type: none"> Zoom in – place two markers at the borders of the graph section you would like to Enlarge. The Zoom icon will change to a magnifying glass with (+) sign . Click the zoom icon. Graph is zoomed-in and the markers disappear automatically. The Zoom icon shows the magnifying glass with the (-) sign . Zoom out – clicking the zoom icon when the icon shows the magnifying glass with (-)  will return the graph to full-view.
	Mathematical Functions – Datahub allows you to apply two different mathematical functions to your data. This drop down menu will allow you to select either: <p><u>Linear Regression</u> – display the best linear line to fit the graph between two markers. A pop up box will display the mathematic formula for the line between the two markers in the $Y = aX + b$ format.</p> <p><u>Quadratic Regression</u> – display the best parabolic line to fit the graph between the two markers. The pop up box will display the mathematic formula for the line between the two markers in the $Y = aX^2 + bX + c$ format.</p>

How To...

	To set a marker on your graph line, simply move the mouse pointer on the line graph, where you would like to place a marker, and double click the mouse left key.
	Activities – open pdf files of lesson plans.

Other Software Features

Marker –set markers on your graph by placing the mouse pointer on the line graph and double clicking mouse left key. In addition to the marker, a pop up bubble with the specific value at that point is added. You can move the marker along the line by pressing the mouse left key and sliding the marker along the line. The values within the bubble will change as you move the marker. You can also relocate the bubble by clicking and dragging it anywhere within 360 degrees of the marker. To remove a marker, place the mouse pointer within the marker bubble, right click the mouse and select delete.

Annotate – add labels to your graph to explain what is happening on the graph by mouse double click anywhere on the graph window **except** on the data line. A pop up bubble is placed on the screen. To add text (or to delete an annotation), double click the mouse when the mouse pointer is within the bubble area. A dialog is opened and you can type in text. Click **Done** and the text will appear in the bubble. To move the entire annotation, place the mouse pointer on the bubble tip. Hold the left mouse key and drag the annotation. *At this point in time*, it is not possible to add an image to an annotation using Datahub on a Chromebook laptop.

Display Options – found in the upper right hand corner of the screen and allow you to select the format for data display. These include:









Line graph – if there is data of more than one type displayed, you can choose which lines to view by clicking on the legend in the upper right hand corner. In the pop up box, you can select line color, format (line or individual data points) or choose to hide a particular set of data.

Bar graph –bar graphs will show data from only one sensor at a time. To change which data is displayed, tap on the legend in the upper right hand corner. In the pop up box, you can select which set of data to display or change the color of the data currently displayed.

Table – show data that in a chart format. This can be used to view exact measurements for every collection.

Meters – for this display type – the Datahub must be connected to the Chromebook. There are several different options for meters including;

Full Dial	Half Dial	Vertical Bar	Horizontal Bar	Numeric Readout	Red, Blue & Green Colorimeter Bulb
					

How To...

To change between different meter types or to adjust the range of a meter, click on the meter to open a pop up window where you can select options.


Map (Google map) – this allows you to overlay a Google Earth map of the area where your experiment was conducted with the data from the sensors in that experiment. You must have an Internet connection for this to work properly, and a GPS must be one of the sensors selected for this data collection. Data is displayed as colored dots, with different colors representing different measurements. If multiple sensors were used, right click the legend on the far right to select which sensor to display. Click an individual dot on the map to see a pop up with the exact measurement for that data point.

Using the Datahub Sensors

When using any of the Datahub sensors, be sure to rotate the outer ring so that the sensor is exposed to the environment during the experiment unless otherwise noted.

Accelerometer (Physics)

No additional set up is needed for the Accelerometer. This sensor measures changes in velocity over time. For acceleration measurements to have any meaning, they must be given in reference to a specific coordinate frame. Acceleration is measured in meters/second².

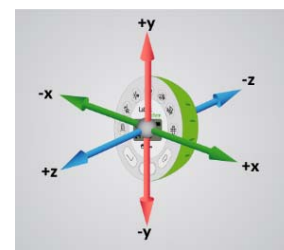

Air Pressure (Gensci, Physics & BioChem)

No additional set up is needed for the Air Pressure sensor, but be sure that the outer ring of the Datahub is rotated to a fully open position to ensure that the sensor is exposed to the environment. Air pressure is measured in kilo Pascals (kPa) which is a metric measure of force per unit area, defined as one newton per square meter. Standard air pressure on Earth is 101.325 kPa.

Ambient Temperature Sensor (All Datahub models)


No additional set up is needed for the Ambient Temperature sensor. This sensor is best when used for data collections longer than 30 minutes. Because it is an internal sensor it may take up to 15 minutes to equalize to the environmental temperature. The Ambient Temperature sensor is primarily used for long term, environmental temperature measurements. The Ambient Temperature sensor can be set to measure in either Celsius or Fahrenheit. To set this, use the configurations menu on the Datahub or your computer.

Barometric Pressure Sensor (Enviro & BioChem)

No additional set up is needed for the Barometer, but be sure that the outer ring of the Datahub is rotated to a fully open position to ensure that the sensor is exposed to the environment. Barometric pressure is measured in millibars (mb). Standard sea level pressure is 1013.23 mb

Colorimeter (BioChem & Enviro)

Before each use calibrate the colorimeter with the solvent you will use during the experiment to help minimize experimental error. To calibrate, fill the cuvette ¾ full with solvent (often distilled water) and place the cuvette in the output. Press and hold the colorimeter button on the Datahub until you hear a long beep. Once calibration is finished remove the cuvette from the colorimeter. To take a data reading, fill the cuvette ¾ full with solution from the sample. Be sure to ALWAYS hold the upper part of the cuvette to avoid leaving fingerprints on the cuvette. Insert the cuvette into the appropriate space and, keeping the top of the cuvette pointed upwards; rotate the outer protective ring shut over the cuvette.

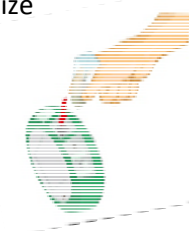

How To...

Conductivity Sensor (BioChem)

Prior to first use, soak the conductivity sensor in distilled water for 30 minutes. Connect the probe to the Datahub and dip in a 0.1 M KCl solution or another standard with a known value. The cell portion of the probe must be fully immersed in the known solution. Stir with the probe to remove any air bubbles trapped within the cell chamber. No air bubbles can be present when measuring. If a 0.1 KCl solution is used to calibrate the meter adjust the meter reading according to the standard value printed on the bottle. After calibration, rinse the probe thoroughly with distilled water.

Submerge the end of the probe in the solution being measured, being sure to completely submerge the cell membrane. Shake the probe well to remove any air bubbles trapped within the cell chamber and check the readout for the solutions conductivity. While the conductivity probe does not tell you which specific ions are present in the sample, it does quickly confirm the total concentration of ions in that sample. Do not touch the probe's cell surface with any hard object. If the cell surface becomes contaminated, soak the probe cell portion in a light detergent and a mild acid for about 15 minutes each and then rinse thoroughly with distilled water. The Conductivity Probe is used to measure the ability of a solution to conduct electrical current. This is done by measuring the concentration of dissolved solids which have been ionized in a polar solution like water. Conductivity is measured in micro-Siemens (mS).

Current Sensor (Gensci & Physics or External Sensor on any Datahub)

Rotate the outer ring of the Datahub to expose the Current Sensor connection port and plug the smaller end of the banana cables into the appropriate ports (red: red, black: black). When using the Current sensor with a Gensci, Enviro or BioChem Datahub, plug the external sensor into the external sensor micro USB port and then plug the smaller end of the banana cables into the appropriate ports on the external sensor. Wait for the Current sensor to auto-calibrate and show a zero reading before beginning data collection. Datahub measures electrical current in Amperes or Amps (A).

Dissolved O2 Sensor (BioChem & Enviro)

Prior to first use, unscrew the membrane module at the bottom of the electrode. Rinse the membrane module with distilled water and blot dry. Completely fill the membrane module with dissolved oxygen electrolyte solution for Galvanic DO Probes. Rinse the internal anode/cathode element with distilled water and blot dry. Screw the membrane module onto the end of the electrode being careful not to over tighten. Rinse with distilled water. Attach the Dissolved O2 sensor to the input bracket by inserting and turning a ¼ turn. Use an aquarium pump to bubble air through 100 ml of water in a 150 ml beaker to completely saturate the water with oxygen, stirring gently with the DO sensor. Keep the sensor at a 20° angle to avoid trapping air bubbles in the sensor. After 20 minutes, hold the Datahub DO2 key down for three seconds until you hear a long chirp. The Datahub should read 8.25 mg/l which is the reading for water saturated with oxygen at room temperature. To use, insert the sensor into the liquid sample and stir slowly and continuously until the value on the Datahub stays constant. Dissolved O2 is measured in milligrams per liter (mg/l)

Distance/Motion Sensor (Gensci & Physics)

The distance sensor is used to measure the distance between or the change in distance between the Datahub and an object. The distance sensor has a viewing angle of +/-15 degrees, so when using it (especially with moving objects) be sure that there are no other objects within that viewing field that might interfere with the distance reading. The distance sensor is located on bottom of the Datahub with a round plastic door covering it. Be sure to fully open the door before beginning data collection with the distance sensor. The distance sensor measures distance in meters.

External Temperature Sensor (All Datahub models)

Plug the External Temperature sensor into the outlet in the Datahub and insert the tip (at least 2cm) into the item/area where temperature is being measured. The External Temperature sensor is suitable for taking environmental temperatures as well as the temperature of liquids and loosely packed solids like soil or sand. The External Temperature sensor can be set to measure in either Celsius or Fahrenheit. To set this, use the configurations menu on the Datahub or on your computer.

How To...

The External Temperature sensor is also available as an external sensor, which allows for the collection of two different temperature readings simultaneously. The external sensor is plugged into the Universal USB port and then the Temperature Probe is plugged into the sensor.

GPS (Global Positioning Satellite) Sensor (Gensci, Enviro & BioChem)

No additional set up is needed to use the GPS sensor. The GPS unit is located behind the housing of the Humidity sensor. It can measure location or speed of movement over a large area. The GPS sensor can also be used to position other collected data onto a Google Earth map, providing accurate information about where the data was collected. The GPS sensor will report the longitude/latitude, speed/course and time/date of the data collected.

Heart Rate Sensor (BioChem or External Sensor on any Datahub)

The ear clip has an IR light source on one side and a light detector on the other side. Blood flowing through the body changes the skin IR transparency, which is measured by this sensor. The ear clip should be placed on the soft section of the ear lobe (an earring hole may interfere with the sensor). You can also attach the sensor clip to the tip of the index finger or the webbing between the thumb and index finger.

If a clear pulse is not observed in the first 5 seconds of measurement, the IR intensity will be increased and checked again for 5 seconds. If a pulse is still not registered, the IR intensity will increase to the maximum intensity. At this point, if there is not pulse detected, it may mean that the clip is connected improperly. For each pulse, the Datahub will produce a short beep.

There are two possible configurations for the Heart Rate sensor.

On the *Biochem Datahub*, the Heart Rate sensor is internal. Plug the ear clip into the port on the Datahub and place the clip onto your ear lobe. The clip will measure both the overall heart rate and the individual waveforms of each heartbeat. After a pulse is detected, the Datahub will count 16 pulses and then start to display the heart rate in beats per minute (bpm).

When using the external heart rate sensor with a Gensci, Enviro or Physics Datahub, plug the external sensor into the external sensor micro USB port and then plug the ear clip into the external sensor. In this format, the clip will measure the individual wave forms of each heartbeat, but will not show a graph line indicating the overall heart rate.

Humidity Sensor (Enviro, BioChem & Gensci)

No additional set up is needed for the Humidity sensor, but be sure that the outer ring of the Datahub is rotated to a fully open position to ensure that the sensor is exposed to the environment. It is used to measure the amount of water vapor in the air expressed as a percentage of the amount needed for saturation at that same temperature. Humidity is measured as a percentage value comparing these two.

Infrared Sensor (Enviro)

No additional set up is needed for the IR sensor, but be sure that the outer ring of the Datahub is rotated to a fully open position to ensure that the sensor is exposed to the environment. An infrared sensor infers temperature from a portion of the thermal radiation emitted by the object being measured. IR is measured as the heat that is radiated off of objects. IR temperatures are measured in Celsius degrees.

Light Sensor (Gensci, BioChem & Physics)

No additional set up is needed for the light sensor, but be sure that the outer ring of the Datahub is rotated to a fully open position to ensure that the sensor is exposed to the environment. Do not point the light sensor directly at the sun.

How To...

This sensor measures the brightness of light in lux. One lux is equal to one lumen per square meter. It is used as a measure of the intensity, as perceived by the human eye, of light that hits or passes through a surface.

Microphone (Gensci & Physics)

There is no additional set up needed to use the Microphone sensor, but be sure that the outer ring of the Datahub is rotated to a fully open position to ensure that the sensor is exposed to the environment. The Microphone measures sound wave forms in voltage and typically records at a very fast rate of 25,000 samples per second.

pH Sensor (Enviro, Gensci & BioChem)

Attach the pH sensor to the input bracket by inserting and turning a ¼ turn. The pH meter will measure the H ion activity in a solution. The pH sensor is factory calibrated and under normal conditions does not require any calibration. If you need to calibrate this sensor; insert the Datahub pH electrode into a 7.0pH buffer and wait for the pH reading to stabilize. Press and hold the pH key for 3 seconds, until you hear a long beep sound indicating that the pH sensor is calibrated and will show a reading between 6.95pH to 7.05pH. To take a pH measurement with the sensor, completely insert the tip into the liquid sample without touching the sides of the container. After each measurement the pH probe should be cleaned with distilled water. Always keep the wash bottle and distilled water close at hand. If you don't have a wash bottle you will need a syringe of at least 10 ml in order to clean the sensor properly. When not in use, the pH sensor should be stored in the buffer solution. The pH sensor measures the number of Hydrogen ions in a solution and is measured on the pH scale of 0 to 14 with readings in the 0 to 6.9 range being considered basic/alkaline and readings in the 7.1 to 14 range being considered acidic. Readings of 7 are considered to be neutral – neither acidic nor alkaline.

Sound Sensor (Enviro)

There is no additional set up needed to use the Sound sensor, but be sure that the outer ring of the Datahub is rotated to a fully open position to ensure that the sensor is exposed to the environment. The Sound sensor measures sound level in decibels. With the Sound sensor, you will typically use 10/second as the sampling rate.

Thermocouple (BioChem)

Rotate open the outer ring to expose the thermocouple outlet. Firmly push the plug of the sensor into the Datahub with the wider prong of the plug to the right. The thermocouple is used to measure temperatures in situations that might not be well suited to a thermometer. Thermocouples have a wider range of temperature measurement; as much as -200°C to +2500°C. Because they are small and have low thermal capacity, they offer a rapid response. Thermocouples are robust and safe, as well as not being prone to self-heating. The thermocouple measures temperature in Celsius degrees.

How To...

Turbidity Sensor (BioChem & Enviro)

Calibrate the turbidity sensor by placing a cuvette full of 1000 NTU turbidity standard into the sensor. Press and hold the turbidity key for 3 seconds until a long beep is heard. To use the sensor, fill the cuvette to a 75% volume with the samples and measure the turbidity, being sure the cuvette is dry outside. Insert the cuvette into the appropriate space and, keeping the top of the cuvette pointed upwards, rotate the outer protective ring shut over the cuvette.

UV Sensor (Enviro)

Be sure that the outer ring of the Datahub is rotated to a fully open position to ensure that the sensor is exposed to the environment. Point the UV sensor directly at the light source from which you wish to measure UV rays. The Ultraviolet sensor measures UV levels with the UVI or Ultraviolet Index. The UV Index is designed as a linear scale, directly proportional to the intensity of UV radiation that causes sunburn. An index of 0 corresponds to zero UV irradiation, while an index of 10 corresponds roughly to midday summer sun with a clear sky. Measurements greater than 11 are quite common in summer at low latitudes, high altitudes, or in areas with above-average ozone layer depletion.


Voltage (Physics or External Sensor on any Datahub)

Rotate the outer ring of the Datahub to expose the Voltage Sensor connection port and plug the smaller end of the banana cables into the appropriate ports (red to red, black to black). When using the Voltage sensor with a Gensci, Enviro or BioChem Datahub, plug the external sensor into the external sensor micro USB port and then plug the smaller end of the banana cables into the appropriate ports on the external sensor. Wait for the Voltage sensor to auto-calibrate and show a zero reading before beginning data collection. Datahub measures Voltage in volts. The volt (V) is the unit of measurement for the electric potential difference between two points.

Available External Sensors

Some of these sensors come built-in with various Datahub models but are available as external sensors for use with other Datahub models.

Dymo Wireless Force & Accelerometer

This sensor allows for wired (via USB) or wireless (via Bluetooth) communication of data relating to force and acceleration. If using Bluetooth communication (recommended for all motion/acceleration activities), you will need to pair the Dymo with your computer the first time you use the sensor. To do this, turn on the Dymo using the on/off button. Press and hold the Bluetooth button for three seconds to enable Bluetooth communication. You will see the blue light flash 3x. Then right click on the Bluetooth icon  in the lower right corner of the Datahub window and click on "Find more Datahubs and sensors." When the "add a device" window pops up, select your Dymo using the last 4 digits from the sticker on the back of the Dymo to identify it. Use 1234 as the pairing code and click "next." Then right click the Bluetooth icon again and select the Dymo that you just paired.

To use the Dymo, press the on/off key. The LED will flash green and the device is on. To start data collection, push the on/off button again (a 3 second hold on the on/off button will turn the Dymo off). When you are finished your data logging, push the on/off button again to stop logging. You can also use the start/stop button in the Datahub software to start/stop your data logging.

If you wish to Zero Force Offset, turn the Dymo off. Remove any weights from the sensor. Disconnect it from the Datahub software. Turn the sensor on again and press both the Bluetooth *and* the On/Off keys simultaneously until the Bluetooth LED is a constant blue. A flashing red light indicates that the battery is low (if unplugged) or charging (if plugged in). Acceleration is measured in meters/second². Force is measured in Newtons (N). A Newton is unit of force needed to move a 1 Kg mass at the rate of 1 m/sec². Since $F=ma$, $1N=1\text{ kg}\cdot\text{m/s}^2$.

How To...

Voltage Sensor

When using an external voltage sensor with a Gensci, Enviro or BioChem Datahub, plug the external sensor into the external sensor micro USB port and then plug the smaller end of the banana cables into the appropriate ports (red/red, black/black). When using the Voltage sensor with a Gensci, Enviro or BioChem Datahub, plug the external sensor into the external sensor micro USB port and then plug the smaller end of the banana cables into the appropriate ports on the external sensor. Wait for the Voltage sensor to auto-calibrate and show a zero reading before beginning data collection. Voltage is measured in volts (V), the unit of measurement for the electric potential difference between two points.

External Temperature Sensor

When using an External Temperature sensor with any Datahub model, plug the external sensor into the external sensor micro USB port and then the Temperature Probe is plugged into the sensor. Since all Datahub models come with an External Temperature sensor built-in, the USB connected External Temperature sensor can be used to measure the temperature of two substances simultaneously. The External Temperature sensor can be set to measure in either Celsius or Fahrenheit. To set this, use the configurations menu on the Datahub or your computer.

CO₂ Sensor

When using a CO₂ sensor with any Datahub model, plug the external sensor into the external sensor micro USB port. The first time the CO₂ is used, you will need to connect the Datahub to its AC/DC adapter and allow the CO₂ sensor to run for 24 hours, to reach calibration. After that, run the CO₂ sensor for at least an hour before starting your experiment.

The power consumption of the CO₂ sensor is rather high, so it is recommended that you have the Datahub plugged into the external power supply when using this sensor. Otherwise Datahub battery will be empty in several hours and it will stop working.

Heart Rate/Pulse

When using an External Heart Rate sensor with a Gensci, Enviro or Physics Datahub, plug the external sensor into the external sensor micro USB port and then plug the ear clip into the external sensor. Place the clip on the fleshy part of the ear lobe, on a small “pinky” finger or on the webbing of skin between the thumb and index finger. With an External Heart Rate sensor, the clip will measure the individual wave forms of each heartbeat, but will not show a graph line indicating the overall heart rate. The pulse waveform will be measured in Volts (V).

Respiration

Prior to using the Respiration sensor with any Datahub model, wash the silicon mouthpiece and dry well before installing it on the Respiration probe. Then connect the Respiration sensor to the Micro USB socket. Turn on the Datahub. If the airflow through the Respiration sensor is within $\pm 10\%$ around zero, the Respiration sensor will auto calibrate and show a zero reading. Have students insert the mouthpiece and begin breathing through the Respiration probe. Respiration is measured in liters per minute (L/min).

Current

When using the External Current sensor with an Enviro or BioChem Datahub, plug the external sensor into the external sensor micro USB port and then plug the smaller end of the banana cables into the appropriate ports (red to red, black to black). When using the Current sensor with a Gensci, Enviro or BioChem Datahub, plug the external sensor into the external sensor micro USB port and then plug the smaller end of the banana cables into the appropriate ports on the external sensor. Wait for the Current sensor to auto calibrate and show a zero reading before beginning data collection. The Datahub measures electrical current in Amperes or Amps (A).

How To...

Magnetic Field

When using the Magnetic Field sensor with any Datahub model, plug the external sensor into the external sensor micro USB port and use the probe's switch to select the Earth's magnetic field range. Use ± 0.2 mT for measuring the Earth's magnetic field or any other very weak field. Use ± 10 mT for measuring the magnetic field near wires carrying current, inside coils and for constant magnetic fields. Magnetic fields are measures in militesla (mT) units.